

KEEPING YOUR DRINKING WATER SAFE

WATER QUALITY MONITORING USING THE H_2S TEST



Tool for Water Quality Monitoring Using The Hydrogen-Sulphide (H₂S) Paper-Strip Test

The Tool for Water Quality Monitoring Using the H₂S Test, is part of the Keeping Your Drinking Water Safe Community Toolkit developed by Live & Learn Environmental Education. The toolkit is designed to be used by Community Trainers, Health Officers, Community Workers, and Facilitators, to raise awareness about the need to keep water clean and promote responsible attitudes, behaviour and actions to ensure safe and lasting drinking water supplies.

Live & Learn Environmental Education is thankful to those who have contributed to the development of the '**Keeping Your Drinking Water Safe Community Toolkit**'. They include:

- The Australian Agency for International Development (AusAID), for funding the 2-year Pacific Drinking Water Safety Plan Programme, which provided financial support for production of this valuable resource.
- The World Health Organization (WHO) and the Pacific Islands Applied Geoscience Commission (SOPAC) for support, guidance and advice during the development of the toolkit.
- The Institute of Applied Science University of the South Pacific, for reviewing the scientific and technical information presented in this toolkit.
- The Government agencies, in-country partners and people of Fiji, Vanuatu, Cook Islands, Tonga, Palau, Samoa, who provided valuable lessons learned.
- The Water & Sewerage Department, Curriculum Development Unit, National Water Quality Monitoring Laboratory and Ministry of Health, Fiji
- Helti Pasifik Komuniti Program, The Foundation of the Peoples of the South Pacific International (FSPI) and affiliates.

The **Keeping Your Drinking Water Safe Community Toolkit** contains:

- An Introductory Guide containing background information and annexes
- Tool for Conducting a Water Audit
- Tool for Conducting Sanitary Surveys
- Tool on Snapshots to Monitoring Water Sources
- Tool For Water Quality Monitoring Using The Hydrogen-Sulphide (H₂S) Paper-Strip Test
- Tool on Water Awareness and Education
- Tool for Water Management Actions
- Comic and Paper-strip test Instruction Flipchart

This document is an output of a regional programme funded by AusAID, for effective management of drinking water supplies in Pacific Island Countries. The views expressed are not necessarily that of AusAID, World Health Organization and/or the Pacific Islands Applied Geoscience Commission.

Illustrations and Graphics by John Robinson

Design and Layout by Live & Learn Environmental Education

Developed by:

Live & Learn Environmental Education

87 Gordon Street, Suva

Private Mail Bag, Suva, FIJI

Ph: +679 331 5868; Fax: +679 330 5868

Email: livelearn@connect.com.fj




Water Monitoring Using the Hydrogen Sulphide Paper-Strip Test

Water Quality Monitoring using the H ₂ S Test.....	2
How Does the H ₂ S Paper Strip Test Work?.....	4
What Can We Use the H ₂ S Test for?.....	5
How Do We Carry Out the H ₂ S Paper Strip Test?.....	6
Filling in the Result Record Sheet.....	10

Annexes

Annex A : Contact List for the H ₂ S Bottles.....	12
Annex B: Water Quality Monitoring Stakeholders.....	13
Annex C: Technical Report on the H ₂ S Test.....	16

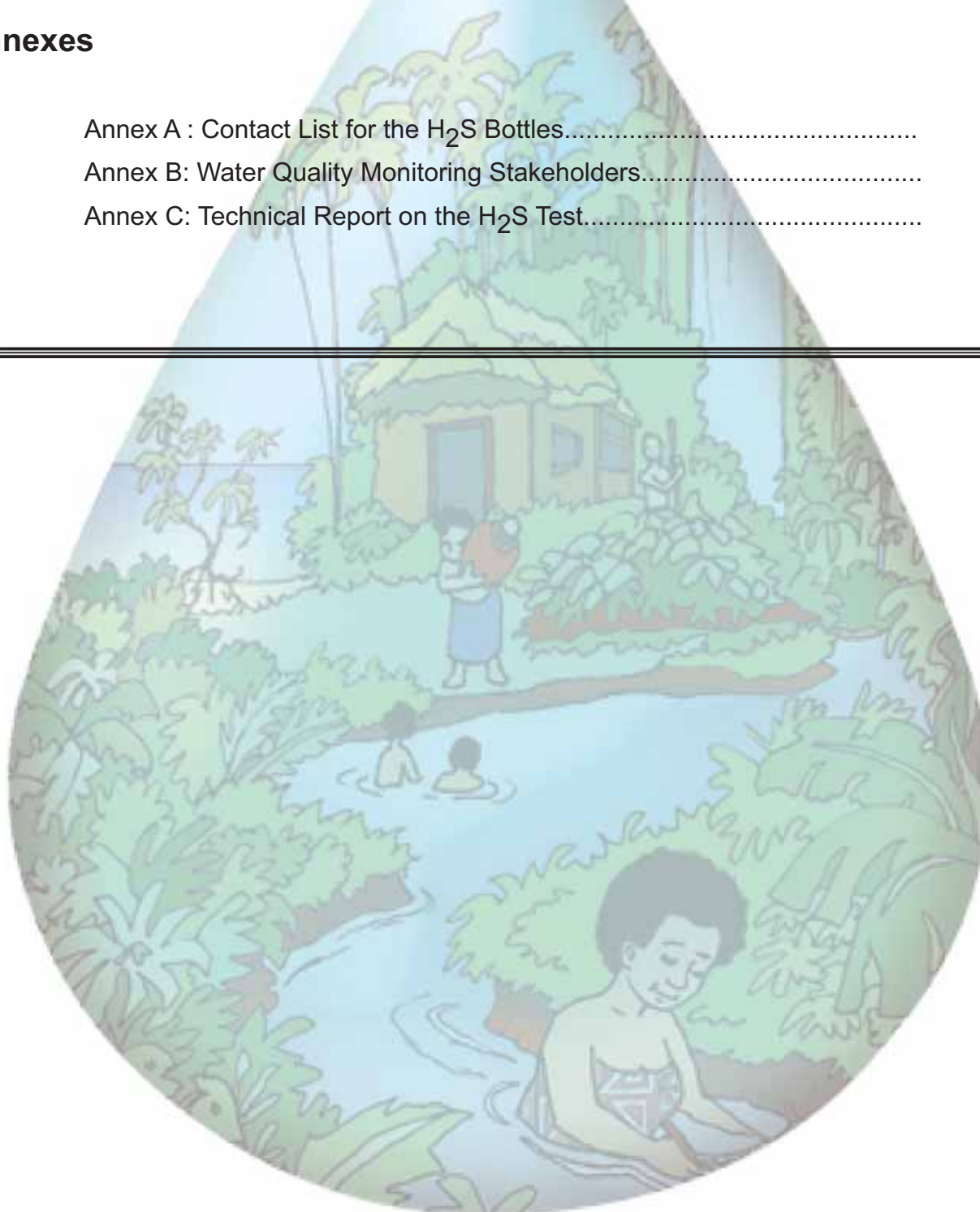


Introduction

Water testing can alert a community to contaminants in time to prevent health problems. Conventional methods of testing require advanced laboratories, highly trained technicians, and expensive supplies - many of which cannot be afforded by most communities or are unavailable in isolated communities.

The Hydrogen Sulphide (H₂S) Paper Strip Test is a drinking water quality test to find out if any of the water being used or stored is already contaminated. It is a very simple test where a water sample is collected, and changes colour if it is contaminated. It's as simple as that!

The Test...

We can test if our drinking water is safe enough to drink by using the **HYDROGEN SULPHIDE (H₂S)** Paperstrip Test

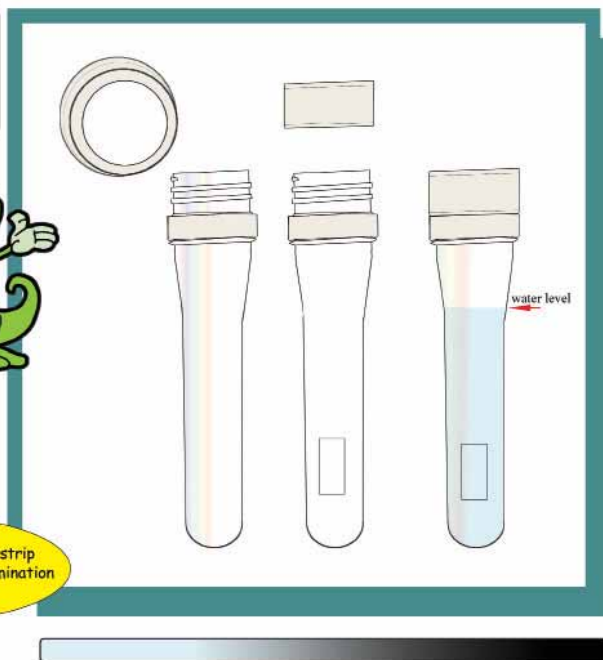

The H₂S Test uses a paperstrip to test for bacterial contamination in our drinking water.




Source: Live & Learn

Getting ready to test one of the wells in Veinuqa village

Testing Our Drinking Water

Contaminated water means that action needs to be taken urgently! The H₂S Paper Strip test can be used regularly to test the quality of water supplies and help people to recognize the risks and take actions if needed.

The H₂S Paper Strip Test was first used in India to test for bacteria contamination in drinking water. Since then many more communities have used it globally and in the Pacific region.

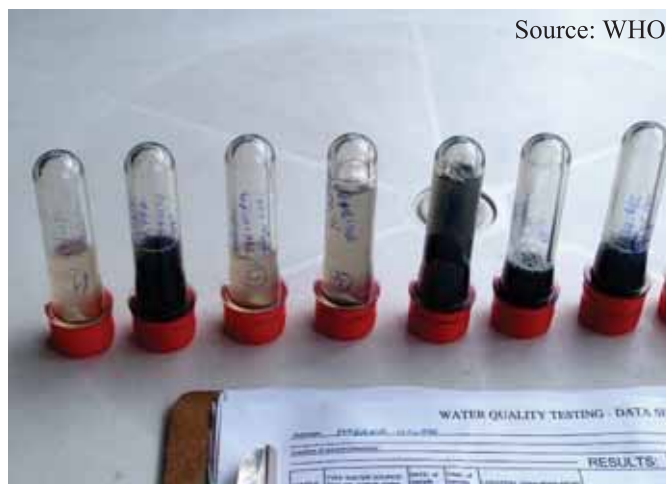

Source: WHO

H₂S Sample bottles and result sheet

The advantages of the H₂S Paper Strip test are, that it is low-cost, does not require samples to be shipped or refrigerated, does not require a laboratory or expensive equipment, and most importantly, it is easy to understand and carry out in the field!

When do we need to test our drinking water?



Whenever we get our water from streams, wells, water tanks or any other source other than a government monitored water treatment plant, the water



needs to be boiled.....



or tested.....



Otherwise it would be the same as taking our water out of our toilet.....



The water may look clean.....

but it is in fact contaminated by dangerous bacteria including the coliform type which can cause the spread of water-related diseases like diarrhoea, typhoid and many more.




How does it work?

The H₂S test is a simple test that will tell us if the water being tested is contaminated within three days (or less). The test identifies if Hydrogen Sulphide (H₂S) is in the sample. H₂S is a gas produced by some harmful bacteria- this is the gas that smells like rotten eggs.

If H₂S is in the water sample, it means that harmful bacteria or viruses from human or animal waste could be in the water.

In order to check for the presence of harmful bacteria in water, a water sample is filled into the test bottle with the paper strip. Chemicals have been mixed into a solution and placed on the paper strip. The paper strip will react with the water sample by turning black if it comes into contact with hydrogen sulphide.

How the H₂S Paper strip Test Works:

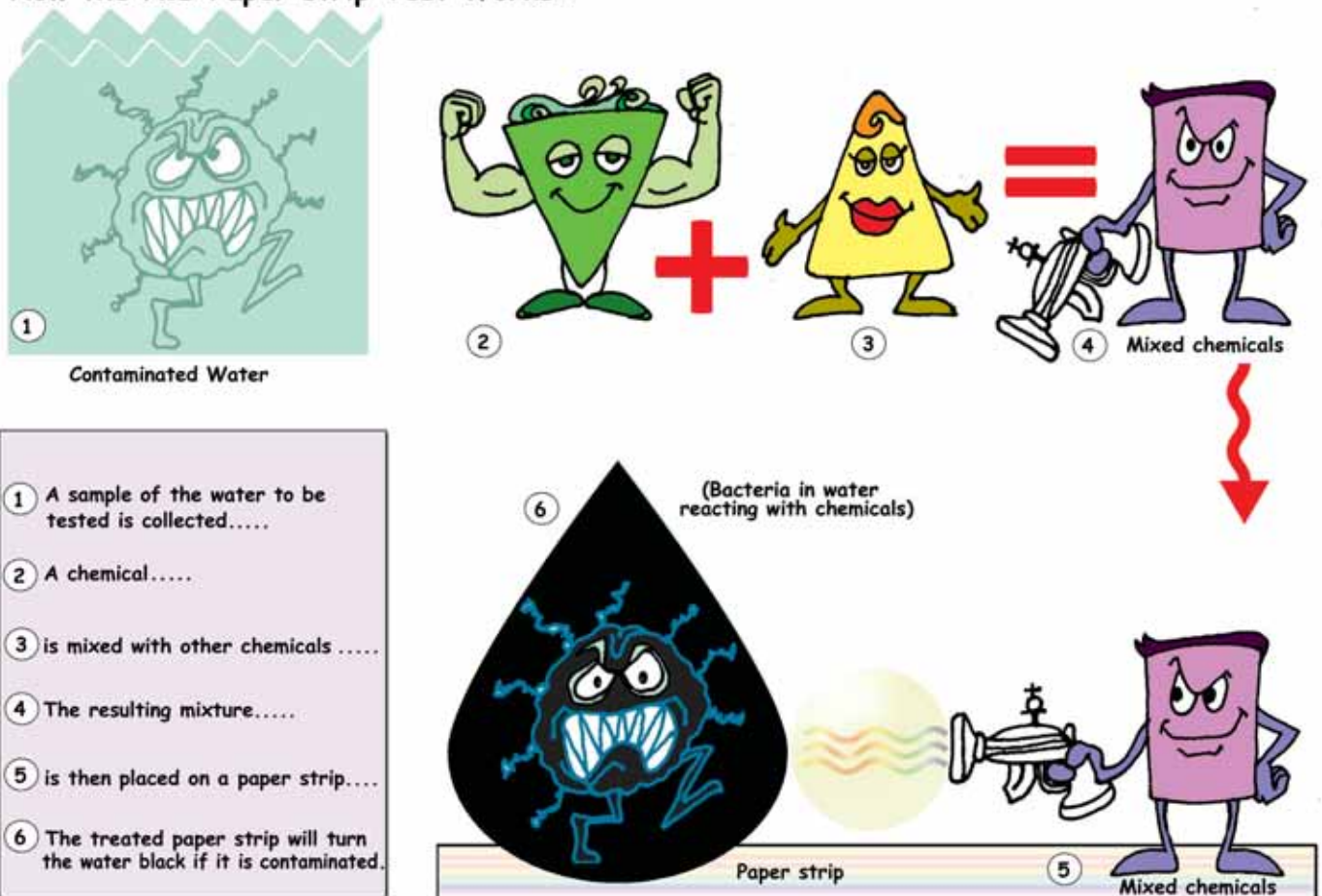

If the water sample or paper-strip turns black, this shows that hydrogen sulphide was produced. This means that it is likely that bacteria are present in the water- that is, the water has been contaminated. This also means that we need to take urgent action.

What can we use the H₂S Test for?

1. For monitoring of rural and outer island water supply systems where it may be difficult to conduct conventional testing due to isolation or a lack of appropriate laboratory facilities.
2. For routine monitoring of reticulated systems; i.e. water that is distributed through a piped system.
3. To identify if there is a need for further analysis of the water sample.
4. To determine the cleanliness of water storage tanks, rainwater cisterns and other household storage containers.
5. To identify sources of contamination or the point in a piped system where bacteria may be entering the water source.
6. To select which spring is best to develop.
7. To check how effective you have been in disinfecting a water source, or to verify that a well has been properly protected.
8. As a tool in health and hygiene education to show villagers how water becomes contaminated and what they can do about it.
9. For monitoring during emergencies and disasters such as cyclones when water-borne diseases are more likely to occur and conventional testing is difficult.
10. To demonstrate how easily hands become contaminated and how easily they can contaminate food and water.



Source: Pacific Islands Applied Geoscience Commission (SOPAC)

The H₂S test can identify whether the water in your storage containers are 'safe' to drink.




Source: Pacific Islands Applied Geoscience Commission (SOPAC)

For example, the H₂S test can be used to demonstrate the effectiveness of washing hand with soap; i.e. to illustrate how bacteria can get from the hands to the mouth and into the body. This can be done by pouring clean water over unwashed hands and testing it, and having others wash their hands with soap and repeating the exercise.

Instructions

Step 1: Filling in the details

- Fill in sample number and date on the round sticker or sticker strip label and stick on the sample bottle (be careful not to get the sticker wet).
- Record your sample number, date, time, location and description of the water source sampled on the Result Record Sheet.
- Record any other information e.g. turbidity, smell, source of pollution, faulty pump etc.

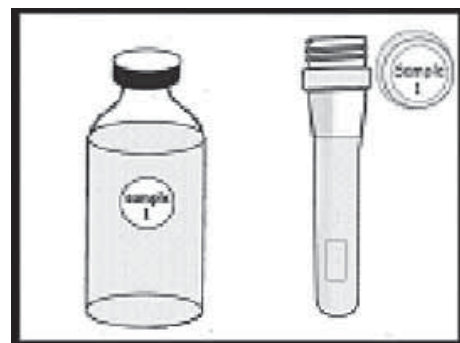

SAMPLE INFORM			
Sample Number	Date	Time	Location- place where sample is
1			
2			


An illustration of a hand holding a pen and writing on a white label strip.

Note!

- Do not open the test bottle until you are ready to fill them with your water sample.
- Make sure that no contamination occurs e.g. by touching the mouth of the bottle.
- Do not hold the cap from the inside.

Step 2: Collecting the Control (Optional)

- Collect a sample of uncontaminated water e.g. distilled water, boiled water, bottled water, water treated with chlorine. This is to be used as the control.
- There may be a slight change in the colour of the sample to a pale yellow or light brown due to the colour change of the reagent. This is normal.

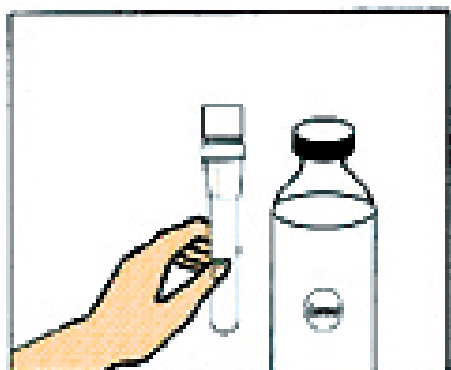

A control is used to compare the colour change in the test samples, and to ensure that the sample bottles are not contaminated before use.

Instructions

Step 3: Collecting the water sample:

A. From the tap





- First clean the mouth or the outlet of the tap with a clean cloth.
- Turn on the tap and allow the water to flow for 15 to 20 seconds.
- Collect sample water from the tap by filling the sample bottle up to the mark. (Fill 10ml or slightly more.)
- Fill the test bottle carefully, this is because the test bottle will fill very quickly to the marked line and may overflow.
- If you do overfill the bottle, do not spill the water out and do not worry. Your result will still be valid.
- Immediately close the sample bottle.



Step 3: Collecting the water sample:

B. From storage containers such as water tanks, and wells or rivers

- Rinse the container to collect the water several times.
- Collect a sample of water from the container by filling the sample bottle up to the mark.
- Close the sample bottle. Make sure that no contamination occurs.

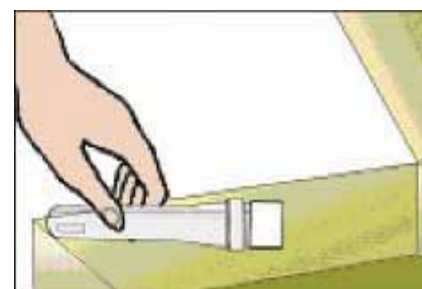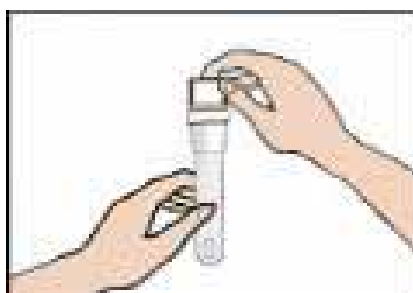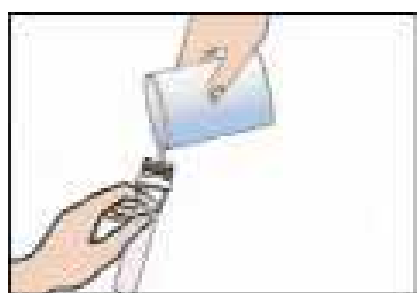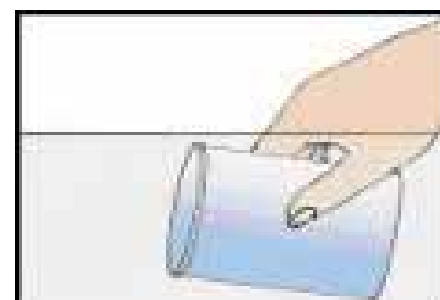


Place all the test samples in a dark place at room temperature.

Do not expose your bottles to direct sunlight. Store in a dark place. The sun's rays can kill the bacteria inside the test bottles and you will not get a true result.

Instructions

Step 4: Checking your results

- Check your test sample at the same time each day for 3 days for changes in colour.
- Use the Result Record Sheet to record the date and time for each observation of your test sample, and your result for each day.
- Compare the colour change of your test sample with that of the control.
- Use the H₂S Colour Code to indicate the degree of contamination.

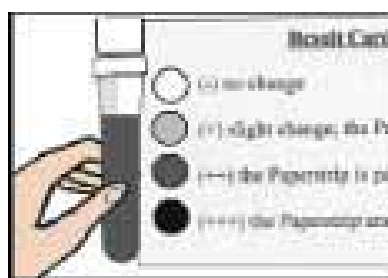


It is important to check your test sample everyday for 3 days, and to keep a good record of your results.


Note!

- Keep the test bottles stored away from children! Do not put it in a place where a child can reach it!
- Do not open used bottles with the water sample. Return the used bottles to get replacements.


H₂S Colour Code

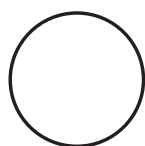

(-) No change

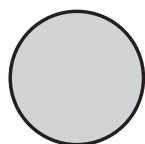

(+) Slight change, the paper strip or water sample has turned grey.

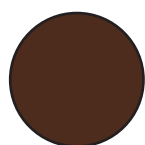

(++) The paper strip or water sample is partially black.







(+++) The paper strip and the water sample are noticeably black.

Instructions

Step 5: Analysing your results. What do your results mean?

- (-) If there is no colour change, this indicates that the water is clean and likely to be free from bacterial contamination.
- (+) If the water sample or paper strip has turned grey, there is a possibility that bacteria is present in the water. **Wait for a few days and check again.**
- (++) If the water sample or paper strip has turned partially black then there is some amount of bacterial contamination in the drinking water. **Conduct a sanitary survey to check your water source.**
- (+++) If the paper strip and the water sample are noticeably black then there is a very high risk of bacterial contamination in the drinking water, therefore, it is not safe for drinking. **Take immediate action! Your water is contaminated!**
- (++++) If there is a fast reaction- that is, the water sample and paper strip turns black overnight, which means that there is a high probability of bacteria present! **Take immediate action! Your water is contaminated!**

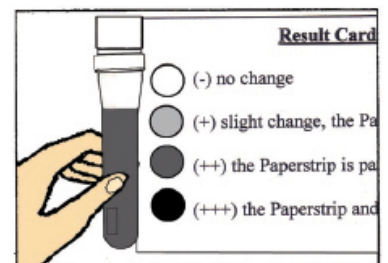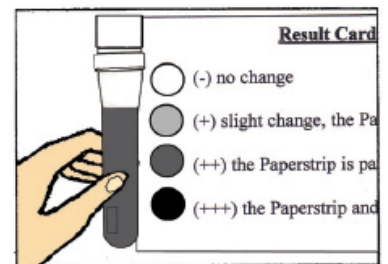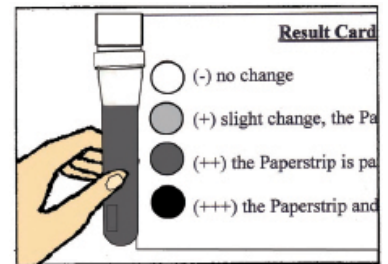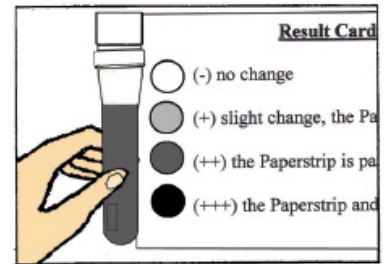

Source: Live & Learn

If your water is contaminated, you should clean out your water storage containers, tanks or well and boil the water before you drink it! Conduct a sanitary survey and check for the source of contamination. Sample the water in your well, tanks and containers again after this to check if you have eliminated the contamination!

Filling in the Result Record Sheet

Every time a Water Monitor is going out for water monitoring, he or she needs to fill in the provided Result Record Sheet. All the relevant details need to be filled.

1. Fill in the address or where you doing the water sampling e.g. Nailega Village, Tailevu.
2. Write in your sample number in the first column.
3. Fill in the type of water source that you are sampling e.g. rainwater tank.
4. Record the date and time of sampling.
5. Identify the location of your sample e.g. the Heilala Community main water tank.
6. In the “Remarks” column, fill in information like the color of the water, the smell, or if there is faulty tap or pipe.
7. Use the colour code to find out your results- e.g. “+” or “++” and record this in the “Results” column. Fill in your observation each day for three days and record the date and time of observation.
8. The “Notes” column can be used for other information like the source of contamination or if there is a latrine/ toilet built within a short distance from the drinking water source.
9. Copy the Result Record Sheet provided on the next page to use and distribute.

Example of the Result Record Sheet

H ₂ S PAPERSTRIP TEST – RESULT SHEET.								
Community Address: _____								
Name of Water Monitor: _____								
RESULTS								
Sample Number	Type of water source: (deep well – borehole, river, rainwater etc.)	Date:	Time:	Location: place where the sample is collected	Remarks	Day 1 Date: Time:	Day2 Date: Time	Day 3 Date: Time:
1								
2								
3								
4								
5								
6								
7								
8								
9								
10								
Notes:								

Result Record Sheet to copy

RESULT RECORD SHEET FOR THE H₂S PAPER STRIP TEST

Community & Address: _____

Name of Water Monitor: _____

Sample Number	Type of water source (deep well, dam, borehole, river, rainwater etc.)	Date	Time	Location (place where the sample is collected)	Remarks (is the water muddy, coloured, contain solids or materials in suspension also an problems at sampling site like a leaking tap, unclean, drainage problems etc.)	Results each day over 3 days (-clear, +grey colour, ++part black, +++very black)
1						
2						
3						
4						
5						
6						
7						
8						
9						
10						

Notes (Distance between water source and possible pollution like a compost pit, septic tank, toilet or farm):

Annex A: Contact List for H₂S Test Bottles

Fiji

Producers:

IAS (USP)

Contact: Prof. Bill Aalbersberg ; aalbersberg@usp.ac.fj

Ms. Arun Pande ; pande_a@usp.ac.fj

Distributors:

WHO (SP)

Contact: Mr. Steve Iddings; IddingsS@wpro.who.int

Ms. Tema Vakaotia; VakaotiaT@wpro.who.int

Cook Islands

Producers:

Ministry of Health

Contact: Ms. Peia Ben; p.ben@health.gov.ck

Department of Water Works

Contact: Mr. Adrian Teotahi; hydro@mow.gov.ck

Distributors:

National Environment Service

Contact: Ms Deyna Marsh; deyna@environment.org.ck

Vanuatu

Producers & Distributors: Ministry of Health

Contact: Ms. Nelly Muru Ham; nham@vanuatu.gov.vu

Tonga

Producers and Distributors: Ministry of Health

Contact: Ms Telesia Talia'uli; ttaliauli@health.gov.to

Palau

Producers and Distributors: Division of Environmental Health (MoH)

Contact: Ms. Joanne Maireng; sengk@palaunet.com

For all other countries, requests for test kits and materials/resources can be sent directly to:

Mr. Steve Iddings or Ms. Tema Vakaotia at World Health Organization (SP) Office, Level 4, Plaza 1, Provident Plaza, 33 Ellery Street, Suva, Fiji. Phone: (679) 323 4100

Annex B: Water Quality Monitoring Stakeholders

COOK ISLANDS

Mr. Adrian Teotahi
Water Quality/GIS Technician
Ministry of Works
Email: hydro@mow.gov.ck
ph: (682) 20034 fax: (682) 21134

Ms Peia Ben
Laboratory Analyst
Health Department
Ministry of Health
Email: p.ben@health.gov.ck
ph: (682) 22664 fax: (682) 22670

FIJI

Mr. Uraia Rabuatoka
Quality Manager
Ministry of Health
Email: uddy_rabuatoka@yahoo.com
ph: (679) 3320066 fax: (679) 3323276

Mr. Nemani Talemaitoga
Technical Officer
Ministry of Health
Email: nemtale@yahoo.com
ph: (679) 3320066 fax: (679) 3323276

Mr. Shelvin Achint Prasad
Technical Assistant
National Water Quality Laboratory, Kinoya
Email: shelvin_achint@yahoo.com
ph: (679) 3392133 fax: (679) 3395495

Mr. Jone Ravulo Tabui
Technical Assistant
National Water Quality Laboratory, Kinoya
ph: (679) 3392133 fax: (679) 3395495

Mr. Sher Singh
Senior Scientific Officer
National Water Quality Laboratory, Kinoya
Email: shersingh@connect.com.fj
ph: (679) 3392133 fax: (679) 3395495

Ms Ranjeeta Ranjani Lata
Scientific Officer
Mineral Resources Department Laboratory
Email: ranjeeta@mrd.gov.fj
ph: (679) 3389441 fax: (679) 3370039

KIRIBATI

Mr. Tianuare Taeuea
Chief Health Inspector
Ministry of Health and Medical Services
Email: btiim@yahoo.com
ph: (686) 28100 fax: (686) 28152

Mr. Tietaake Beia
Water Production Foreman
Public Utilities Board
Email: c/ Itienang Timona itimona6@yahoo.com
ph: (686) 26292 fax: (686) 26106

NIUE

Mr. John Hetutu
Chief Environmental Health Officer
Health Department
Email: environ.health@mail.gov.nu
ph: (683) 4100 fax: (683) 4265

PAPUA NEW GUINEA

Mr. Pius Palma
Principal Officer – Water Quality
PNG Waterboard
Email: ppalma@pngwater.com.pg
ph: (675) 3235700 fax: (675) 3256298

Mr. Raka Jr Taviri
Executive Manager Operations
PNG Waterboard
Email: rtaviri@pngwater.com.pg
ph: (675) 3235700 fax: (675) 3256298

Ms Dianne Kave
Environmental Health Officer
Department of Health
ph: (675) 3013704 fax: (675) 3013604

SAMOA

Mr. Afaesefau Meredith
Laboratory Technician
Samoa Water Authority
Email: afa@swa.gov.ws
ph: (685) 21267 fax: (685) 21298

Mr. Paulo Pemitia Seuseu
Senior Water Quality Officer
Ministry of Health
Email: PauloP@health.gov.ws
ph: (685) 21212 fax: (685) 21106

SOLOMON ISLANDS

Mr. Wycliff Mamou Maebule
Quality Control Officer
Solomon Islands Water Authority
Email: c/ jhoutarau@siwa.com.sb
ph: (677) 23985 fax: (677) 20723

Mr. Jay Semah Kabei
Chief Laboratory Scientist
Ministry of Health
Email: jkabei@simtri.gov.sb
ph: (677) 38871/25513 fax: (677) 25513

TONGA

Mr. Timote Fakatava
Head Chemist
Tonga Water Board
Email: timotefakatava@yahoo.com.au
ph: (676) 23299 fax: (676) 23518

Ms Telesia Talia'uli
Medical Scientist
Ministry of Health
Email: ttaliauli@health.gov.to
ph: (676) 23200 fax: (676) 24210

TUVALU

Mr. Pisi Seleganiu
Water and Sewage Supervisor
Public Works Division
Email: seleganiu@yahoo.co.nz
ph: (688) 20306/20556

Mr. Falealili Feagai
Senior Environmental Health Officer
Ministry of Health
Email: lfaavae@yahoo.com
ph: (688) 20480 fax: (688) 20481

VANUATU

Mr. Erickson Sammy
Water Resources Manager
Department of Geology, Mines and Water Resources
Email: amapelao@yahoo.com
ph: (678) 22423 fax: (678) 22213

Mrs Nellie Muru Wouloseje
Environnemental Health Officer
Ministry of Health
Email : nham@vanuatu.gov.vu
ph: (678) 22512 fax: (678) 25438

RESOURCE PERSONNEL

World Health Organization
Mr. Steven Iddings
Environmental Engineer
Email: IddingsS@wpro.who.int

Mr. Mitesh Mudaliar
Programme Officer – Pacific Water Safety Plans/
Water Quality Monitoring
Email: mudaliarm@wpro.who.int

Guam Environmental Protection Agency
Mr. Angel Marquez
Safe Drinking Water Programme Director
Email: Angel.Marquez@guamepa.net

Mr. Rodolfo Paulino
Chemist
Email: Rodolfo.Paulino@guamepa.net

Institute of Environmental Science and Research
Limited (ESR)
Mr. Andrew Ball
Microbiologist
Email: Andrew.Ball@esr.cri.nz

IDEXX Laboratories Australia
Mr. Brendan Gee
Account Manager
Environmental Products
Email: brendan-gee@idexx.com

Institute of Applied Sciences, the University of the
South Pacific
Professor Bill Aalbersberg
Director
Email: Aalbersberg@usp.ac.fj

Ms Apiame Cegumalua
Fellow – Administration
Email: cegumalua_a@usp.ac.fj

Mr. Usaia Dolodolotawake
Acting Lab Manager
Email: dolodolotawa@usp.ac.fj

Mrs Mereoni Gonelevu
Quality Assurance Officer
Email: Gonelevu_m@usp.ac.fj

Mrs Arun Pande
Senior Technician
Email: pande_a@usp.ac.fj

Mr. Girish Lakhan
Assistant Project Manager
Email: lakhan_g@usp.ac.fj

Ms Ashika Dayal
Graduate Assistant
Email: ashik_d@yahoo.com

Pacific Islands Applied Geoscience Commission
(SOPAC)
Mr. Tasleem Hasan
Water Quality Monitoring Programme Officer
Email: Tasleem@sopac.org

Mrs Arieta N Sokota
Water Project Officer
Email: arieta@sopac.org

Mr. Marc Overmars
Water Adviser
Email: marc@sopac.org

Mr. Kamal Khatri
Water, Sanitation and Hygiene Officer
Email: kamal@sopac.org

Mr. Davendra Nath
Water Safety Plans Officer
Email: davendra@sopac.org

Ms Sanjeshni Naidu
Water Partnership Officer
Email: Sanjeshni@sopac.org

The Hydrogen Sulphide (H₂S) Paper-Strip Test

A simple test for monitoring drinking water quality in the Pacific Islands

Luke M. Mosley¹ and Donald S. Sharp^{2,}*

¹South Pacific Applied Geoscience Commission (SOPAC)

²World Health Organization (WHO), Suva, Fiji Islands*

SOPAC Technical Publication 373

March 2004

**Disclaimer: This document reflects the views of the authors and not necessarily those
of WHO*

Introduction

Many people living in the Pacific Island region are reliant on drinking water from shallow ground water lenses and streams, or from roof catchment systems. These water resources are often scarce and vulnerable to contamination from poorly installed sanitation facilities (Falkland 1999). The majority of the Pacific's population therefore is at risk to waterborne diseases (Prüss et al., 2002). Many islands are too small in size and under resourced to support conventional water treatment facilities and distribution networks. In urban areas, distribution and treatment systems are reasonably common although maintaining the systems in good working order is sometimes difficult. People living in rural areas and outer island groups are most at risk of contracting waterborne diseases, as they do not typically have access to treated water.

Microbiological and chemical testing of drinking water quality should be performed to indicate whether water is safe to drink. Unfortunately, in many Pacific Islands the infrastructure needed to adequately monitor water quality is either non-existent or inadequate. A lack of monitoring is particularly apparent for outer islands and rural areas.

Typically, the key water quality parameter that indicates safety is the absence faecal coliform bacteria in the sample. These bacteria indicate the probability of pathogens (e.g. typhoid and cholera) being present. Sophisticated and costly equipment is required to test for these organisms; i.e. an incubator, filtration apparatus, and chemical reagents, which must be stored under refrigeration. The cost of one test can be more than NZ\$30, depending upon the type of test and method used. In addition, samples for coliform analyses must be kept chilled and delivered to the laboratory within 6 h. in order to adequately preserve the sample. In all cases, the time elapsed between collection and examination should not exceed 24 h (APHA, 1995). Although there are several commercially available portable kits that make it possible to carry out on-site water quality testing, these are usually costly and require technical expertise to operate (Bartram & Balance 1996).

An alternative low-cost test for faecal contamination in drinking water which is simple to use and easy to interpret is the hydrogen sulphide (H_2S) paper-strip test (Manja et al. 1982).

Aim of this Report

The aim of this report is to provide information on the scientific basis, manufacture and use of the H_2S paper strip test in the Pacific Islands. Suggestions are given on how the test could be utilised for rural, outer island and community-based water quality monitoring.

Indicators of faecal pollution

Untreated or improperly treated drinking water may contain microorganisms of faecal origin that are pathogenic (disease causing) such as those that cause cholera and typhoid fever. The presence of pathogens in drinking water is usually due to human and animal waste entering the water source. The sanitation facilities that are used predominantly in rural/outer islands of the Pacific are septic tanks and pit latrines, which do not provide sufficient treatment to remove pathogens. The waste outflow from these types of facilities, in certain soil conditions can travel several hundred metres underground (Dillion 1997). Animals (e.g. pigs and cows) in the area of an unprotected water supplies can also cause serious contamination and pose risks to public health.

It is difficult and expensive to test for the pathogenic organisms that may be present in contaminated drinking water. Therefore indicator organisms are used to determine the risk that these organisms might be present in drinking water. Indicator organisms are always present in high numbers in faecal material, whether or not pathogenic organisms may be present. A high level of indicator organisms in a water sample indicates a high risk that pathogenic organisms might also be present. The most common indicator organisms used to determine bacteriological water quality are total and faecal coliforms.

The coliform group of bacteria, along with many other naturally occurring bacteria, inhabit the intestinal tract of warm-blooded animals, including humans, and are discharged in their faeces. Faecal coliform presence indicates that water is contaminated with faecal matter and is not safe for drinking purposes. In the tropics, coliforms are not an ideal indicator as they can occur naturally and reproduce in soil and water at the ambient temperatures (WHO 1996). Other indicator organisms are sometimes used which are in the Enterococcus bacteria group such as faecal streptococci (WHO, 1996), and Clostridium perfringens. The problems noted above with the sophisticated testing procedures and equipment required for the analysis of the above indicator organisms make their use difficult in rural areas and on outer islands.

Another less commonly used indicator is sulphide-reducing bacteria and the following sections outline the use of a low cost test for these bacteria in drinking water called the hydrogen sulphide (H₂S) paper-strip test. There are many advantages of this test for use in rural and remote Pacific Island communities particularly where conventional monitoring is not possible or too expensive.

Background and scientific basis of the H₂S test

In 1975, Allen and Geldreich showed that the presence of coliforms in water was also associated with hydrogen sulphide (H₂S) producing organisms. In 1982, Manja et al. developed a simple paper-strip method to screen for bacteriological contamination of potable waters. This study and several subsequent studies, have found that the H₂S test gave generally good agreement with the standard Most Probable Number (MPN) and membrane filtration methods commonly used for determining the presence and number of coliform and faecal coliform organisms (Hazbun & Parker 1983; Dutka 1990; Castillo et al. 1994; Martins et al. 1997; WHO 2002). As noted in a multi-country intercomparison study summarised by Dutka (1990), this test is “an ideal tool for testing rural and isolated drinking water supplies”.

Bacteria can produce hydrogen sulphide through the anaerobic catabolism of cysteine, an amino acid containing the sulfhydryl group, or by the use of elemental sulphur or some oxidised sulphur compounds as the terminal electron acceptor in their metabolic processes. All members of the Enterobacteriaceae group are capable of the former while the latter occurs in dissimilatory sulphate-reducing bacteria. The H₂S test uses a medium with thiosulphate as a sulphur source and ferric ammonium citrate as an “indicator,” only certain enteric bacteria will produce hydrogen sulphide resulting in the development of a black precipitate. Hydrogen sulphide is produced by the reduction of thiosulphate and then reacts with the ferric salt to form an insoluble black ferrous sulphide precipitate. Members of the Enterobacteriaceae group such as Salmonella, Citrobacter, Clostridia, Klebsiella and Proteus are all able to produce hydrogen sulphide in such a medium. Some other non-gut bacteria can reduce thiosulphate into hydrogen sulphide in anaerobic conditions. These bacteria are not typically present in drinking water. The presence of the iron as an indicator in the H₂S medium, would inhibit some naturally occurring bacteria from producing hydrogen sulphide.

The Codex Alimentarius Commission recommends the use of sulphite reducing anaerobes as an indicator for testing bottled natural mineral water (CAC/RCP 48-2001). A report by WHO (2002) did not recommend the use of H₂S bacteria for routine monitoring of water supplies due to the possibility of false positives from naturally occurring sulphite reducing bacteria. Nevertheless, these bacteria can be a valuable tool in that they show a lack of sanitary protection somewhere within the system and indicate a need for further investigation and/or treatment.

Basics of the H₂S paper strip test

The H₂S test is recommended for testing drinking water derived from surface water, boreholes, and rainwater sources for faecal contamination¹.

The reagents used to make the H₂S paper strip test are common laboratory chemicals. By adding a measured amount of “boiled” water and a common liquid detergent to the reagents, a measured amount is impregnated on a piece of absorbent paper and dried in a low-temperature oven. The dried paper strip is placed in a clear small plastic or glass bottle or tube. A water sample is collected in the container containing the reagents and stored in the dark at room temperature for about 3 days. If the sample contains hydrogen sulphide producing organisms, the pad and water turn black in. The black colour and the rotten egg smell of hydrogen sulphide clearly indicate that there is a problem. With such an indicator it is not difficult to convince uneducated villagers that the water may not be safe to drink.

Advantages of H₂S test for use in the Pacific Islands

The advantage of this test over other more sophisticated analyses like the total and faecal coliforms (membrane filtration and Most Probable Number) methods is that it is:

1. Low in cost, the cost of reagents for one test is estimated at NZ\$0.08 (see Appendix 4). All other materials used can be found locally.
2. When making up and using the test it is not necessary to have access to a laboratory or expensive equipment like an autoclave or incubator. Only a simple balance to weigh the media, pipettes, and a method of sterilizing the kits (hot oven, autoclave, UV light) are needed.
3. Does not require samples to be shipped or stored under refrigeration.
4. Samples are incubated at room temperature.
5. Is very easy to use in the field as it consists of only a sample tube.
6. Simple for non-technical people to understand as a clear colour change is observed.

For these reasons the test can be distributed to households/communities so they can test their own water. With sufficient public education on the test, there should be no need to go back and tell them that their water is safe or contaminated as they are conducting the test themselves. If results indicate high risk, households would be instructed to treat their water to make it bacteriologically safe before drinking.

¹The test is not currently recommended for use in testing seawater.

Disadvantages of H₂S test for use in the Pacific Islands

1. Sulphide reducing bacteria (responsible for production of H₂S) are common in the intestinal tracts of most animals making them good indicators for faecal contamination. However, there are some bacteria within the group that occur naturally around thermal vents, vegetation undergoing bacterial decomposition, etc., which may yield a false positive test result (WHO 2002). For most drinking water supplies and all rainwater cistern systems, such 'false positives' would not be expected. As mentioned above there are members of the coliform group that are naturally occurring as well.
2. Some H₂S producing bacteria such as *C. perfringens* are spore formers and hence they may be present long after a pollution episode has occurred (WHO, 1996). However, pathogens may also survive for long periods of time in the tropics (Dillion 1997).
3. Another criticism of the test has been its use as a presence/absence test. The number of indicator organisms in a water sample can indicate the degree of contamination and therefore relative risk to public health. The H₂S test just indicates whether there is a risk, not the degree of risk. However, the speed of the reaction (color change from clear to black) indicates bacterial density. The faster the reaction, the greater the numbers of organisms present. If necessary, estimates of the concentration of bacteria of faecal origin can be made by controlling the volume of the sample used, or by using a three tube or five tube series Most Probable Number method.

Making the H₂S paper strip test

1. Any type of glass bottle or plastic tube with a volume of between 15-200 ml, which has a heat resistant cap/lid, can be used. The bottles or tubes are first cleaned by washing in detergent, rinsing with tap water, soaking in a 5% bleach solution overnight and rinsing with tap water and drying in air or in an oven.
2. If no volume marks are present on the bottles, they can be marked at 10ml, 20ml or 100ml volume, or any volume in-between – depending upon the bottle size. The authors typically use bottles marked at a 10ml volume. This volume calibration is typically done by measuring the required (i.e. 10, 20, 100mL) of water into a graduated cylinder or other measuring device, pouring the measured volume into a sample bottle, standing the bottle upright and then making a mark on the bottle where the water level is. By lining up this bottle with another bottle, the other bottles can be marked in

approximately the same place. A glass marking pencil, permanent ink pen or tape can be used to mark the desired volume.


3. The medium used in the test is prepared from the following chemicals, which are dissolved into distilled or boiled tap water. Shake or stir the mixture to dissolve the chemicals.

H₂S Media Formula

Bacteriological peptone.....	40.0g
Dipotassium hydrogen phosphate.....	3.0g
Ferric ammonium citrate.....	1.5g
Sodium thiosulphate.....	2.0g
Citrate (optional but increases sensitivity).....	g?
Liquid detergent (e.g. Teepol)	2.0ml
Water (distilled or boiled tap).....	100.0ml

4. Taking absorbent paper, filter paper, non-toxic paper toweling, gauze, absorbent pads used for membrane filtration, or any other type of absorbent material², place a measured quantity of media onto the paper. Each paper strip for a 10ml test sample needs to contain 0.5ml of media (50ml sample will use 1ml of media and a 100ml sample will require 2.5ml of media etc³). Large adsorbent pads can be cut to a size that has absorbed 0.5ml of media. For example in the picture below, the pad is ready to be cut into eight paper strips, therefore 4mL of the media will have been poured onto the pad (0.5 mL per strip).

Percentage Distribution of Water on Earth

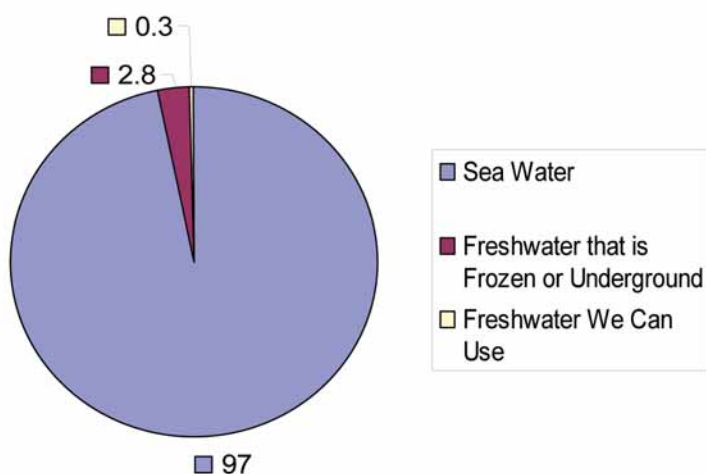

²Coasters used in bars work well (if no black ink is used).

³A pipette graduated at 0.1ml intervals may be necessary for this.

5. The next step is to dry the strips in an oven at about 55°C. A conventional household oven on low temperature can be used. These reagent-impregnated strips can be stored dry (in an envelope or preferably a zip locked bag) for several months – until ready for use.
6. Before conducting the test, a strip or strips are introduced into the appropriate (clean) sample bottle.



7. Bottles should next be loosely capped and sterilized by various possible means:
- Plastic and/or glass tubes can be placed in a hot air oven at about 120°C for 60 minutes.
 - If the tubes are clear plastic they can be sterilized under UV-light for at least 30 minutes.
 - If the tubes are autoclavable-glass (pyrex), sterilization can be done with an autoclave for 30 minutes.
 - Tubes can be placed in a simple pressure cooker for 15 minutes at 115°C.
 - It can also be done by steam (in a rice steamer) for about 30 minutes.

Sterilization in a hot-air oven

Following any of the heat treatments, the tubes or bottles are then allowed to cool and the caps or lids tightly sealed. The tubes or bottles should be stored in a dark place until ready for use. Experience has shown they can be stored for at least 5 years in this manner.

Alternatively, the media can be prepared in pre-weighed, dry form and stored until ready for use. A central laboratory could make the impregnated pads (filter paper) and provide these in sealed plastic “zip-locked” bags or envelopes, or provide ready-made tubes or sample containers. In dry form, these reagent-impregnated pads can be stored indefinitely. The only step required of the community is to insert the reagent coated paper strip in sample containers 15-200 ml in size, and sterilize.



Sampling Procedure:

A basic instruction sheet for distribution to individual households on sampling and interpreting results is shown in Appendix 1. The following procedures are for conducting a survey of water quality using the H₂S test.

1. At the time of sampling, label each container with a sample number.
2. Write the sample number, date and time of collection on the special report form (see Appendix 2). Include on the sheet additional information on the type of water sampled (well, surface, rain water, treated supply, etc.) and exact location of collection, such as “at the tap nearest the borehole.” Under remarks note if the water was visibly turbid or has any other characteristic that should be noted.
3. At the bottom of the report form there is space reserved for “Notes.” Record any observations that may have influenced the quality of the water sampled. For example, distance to a nearby source of pollution, faulty pump, or state of sanitary protection (if a well or spring).
4. Standard procedures specify that if the sample is from a tap, flame the mouth of the tap to eliminate the chance of accidental contamination (a false positive), then, let the water run freely for about 15-20 seconds. Place the opened H₂S sample collection bottle under the tap and collect the appropriate amount (e.g. up to 10mL calibration mark) being careful not to contaminate the cap. It should be noted however that samples should not be collected from taps that are leaking and flaming the tap is not necessary if you are testing the quality of the water as it is actually consumed.
5. If the water to be sampled is from a storage container, tank or cistern; a natural flowing water body like a spring or stream; or from a dug well, use the utensil that is normally used by the consumer or water collector to collect the sample, rinsing it several times before collecting the sample.
6. Every day of sampling, a control sample should be collected and analysed. This is a sample that is known to be uncontaminated, such as boiled water, commercially bottled water, or water treated with chlorine. The control sample is used as a benchmark to compare color change in the test samples and to insure that the sample bottles have been properly sterilized prior to use (Note: There will be slight change in the color of the sample to a pale yellow or light brown due to the color of the reagent, which is normal).

Reading and Interpreting Results:

1. After sampling, place all test samples in a dark place and incubate at room temperature for a total of three days. Every 12-18 hours examine the samples for changes in color. The date and time of each observation is recorded on the report form and the observations are recorded as follows: (-) = no change; (+) = slight change, the paper strip or water has turned gray; (++) = the paper strip is partially black; (+++) = the strip and the water sample itself are noticeably black.



2. As noted above, a color change indicates the presence of bacteria of faecal origin. The speed of the reaction will determine the density of organisms present; i.e. the quicker the reaction the higher the number of faecal organisms presence. This can also be interpreted in terms of a risk factor. For example, a slight color change (+) on day three indicates a lesser risk than a strong (+++) change on day 1.

3. To determine actual risks to health, H₂S test results must be considered in parallel with the results of a sanitary survey. An example of a sanitary survey form for rainwater tanks is shown in Appendix 3. For example, if a drinking water well is unprotected and the results of the H₂S test are positive on the first day, the users should be informed that a risk to health is likely, and steps must be taken to disinfect the water. Usually after seeing the results first hand, the user understands that the water supply in question is not suitable for drinking purposes. In such cases the users are generally receptive to taking corrective action; i.e. to protect the well from contamination, or to learning about disinfection (treating water to kill bacteria).

Various household disinfection techniques could be recommended such as boiling, adding a few drops of chlorine bleach (4 drops per litre), and/or putting the water in a clear plastic bottle and exposing it to full sunlight for a minimum of 4 hours (e.g. 10:00am-2:00pm). Two H₂S tubes could be distributed to each household, with instructions to fill one with untreated water and the other with water that has been treated. After the 2-3 days of incubation, no color change should occur in the treated sample, which clearly shows that the organisms that caused the untreated sample to turn black have been deactivated.

How the H₂S test can be used:

This low-cost test has several useful functions⁴. This test can be used:

1. For monitoring of rural and outer island water supply systems where it may be difficult to conduct conventional testing due to a lack of appropriate laboratory facilities. This, along with sanitary survey data (information identifying conditions that may lead to the contamination of a water supply source) would make it possible for communities to monitor their own water supplies without having to rely on central laboratory services. If community-based programmes of this nature can be established, outer islands could carry out their own water quality monitoring and surveillance programmes and initiate corrective action when needed. This would help to protect public health in these areas and result in a substantial cost savings to island governments.
2. For routine monitoring of reticulated systems; i.e. water that is distributed through a piped system. If a positive result is observed, another sample can be collected for further analysis by conventional means e.g. for faecal coliform enumeration;
3. To determine the cleanliness of water storage tanks, rainwater cisterns and other household storage containers;
4. To identify sources of contamination or the point in a reticulated system where bacteriological contamination is being introduced.
5. To select which spring is best to develop;
6. To determine effectiveness of disinfecting a water source, or to verify that a well has been properly protected;
7. As a tool in health and hygiene education to show villagers how water becomes contaminated and what they can do about it. Communities would also have evidence to alert the relevant authorities that water which is supposedly treated is still contaminated.

⁴Keep in mind that this is a tool to illustrate bacteriological contamination and is not a standard method that is admissible in legal proceedings. The major usefulness of this test is its application as a low-cost educational tool.

8. For monitoring during emergencies and disasters such as cyclones where conventional testing is difficult. For example, following a cyclone, thousands of kits could be locally manufactured and distributed in a short time period (few days) to individual households by community health workers along with printed material on their use. As most water sources on outer islands are localized small community or individual sources, time and resources could be focussed on distribution of the kits to the maximum number of people. This would mean that many more drinking water supplies are tested compared to if conventional methods were used by a water or health agency. This should enable better protection of human health following disasters.
9. To identify sources of pollution entering streams and rivers by comparing differences in incubation times of samples collected at regular intervals along the stream or river;
10. To demonstrate how easily hands become contaminated and how easily they can contaminate food and water. For example, it can be used to demonstrate the effectiveness of washing hand with soap; i.e. to illustrate the faecal oral route of disease transmission. This is done by pouring clean (boiled and cooled) water over unwashed hands and testing it, and having others wash their hands with soap and repeating the exercise; and,
11. To determine if a food contact surface is contaminated with an a common foodborne pathogen like Salmonella by swabbing the surface with a sterile swab and inserting the swab in an H₂S tube containing sterile water.

Commercial availability of H₂S test

Although the test is very simple to make and this should be possible in most countries, the test may also be purchased when manufacture is not possible.

1. H₂S tests or chemical reagents may also be able to be obtained on a 'cost of manufacture only' basis from the

Institute of Applied Sciences,
University of the South Pacific,
Box 1168,
Suva,
Fiji Islands.

pH: (679) 3212967 or 3212965

Fax: (679) 3300373

See website for email contact details: www.usp.ac.fj/ias

2. HACH chemical company make a H₂S test called the Pathoscreen test.

Validation of H₂S test for use in the Pacific Islands

The H₂S test has been used in several Pacific countries but no detailed comparisons have been published where this test method is compared against other methods of assessing the microbiological quality of water. For this reason we undertook laboratory and field testing and validation procedures to ensure the test was suitable for use in the Pacific Islands.

Laboratory Testing

In order to determine how this test can be developed for use in developing countries, a series of experiments were conducted on three naturally occurring water types which are commonly used as drinking water sources:

1. Water from a large river (currently used as a water supply) flowing through a rural catchment.
2. Water from a small creek flowing through a semi-urban area.
3. Water from a typical household rainwater cistern system found in Fiji.

The intent of these experiments were to determine how well the test correlates with other traditional water quality tests and whether or not the time it takes for a reaction to occur correlates with bacterial density and relative risk. All analyses were undertaken using validated methods in the microbiological laboratory at the Institute of Applied Sciences, University of the South Pacific, Suva, Fiji.

The table below shows the time taken for H₂S development compared to bacteria counts (Colony Forming Units per 100 mL, CFU/100mL) using conventional methods for total (TC) and faecal (FC) coliforms. The time taken for H₂S development is separated into the time taken for initial (grey colouration, +) and final (dark black colouration, +++) colour development. The creek water was the most contaminated water with very high counts of total and faecal coliforms, and this water also took the shortest time to turn black. The rain water took longer (92 h) to turn fully black and had low levels of faecal contamination. We also examined the use of the H₂S method to determine bacteria counts using a Most Probable Number (MPN) method (serial dilutions made of sample and these dilution placed in separate H₂S tubes). The results agreed well with the trends in the other results indicating that, if necessary, the H₂S test could be used to estimate bacteria numbers.

Site	Time for H ₂ S development	Time for H ₂ STC development	FC	MPN H ₂ S
	initial colour	full colour	CFU/100 mL	CFU/100 mL
River water	40 hours	59 hours	480000	62
Creek water	23 hours	25 hours	3820000	2700000
Rainwater	42 hours	92 hours	490000	1

The same samples were tested for faecal streptococci (FS, Enterococci), Salmonella (2 different types of test) and Clostridium perfringens. The faecal streptococci results were similar to the faecal coliforms which is not unexpected as both these indicator bacteria groups would be expected to be found together in faecally contaminated water. *C. perfringens*, a definite indicator of faecal pollution, was also found in the river and creek samples but not the rain water. Salmonella results were variable with one test showing positive for creek water and one test showing positive for river water. Salmonella would be expected to be at much lower levels than the other types of bacteria so the statistical probability of finding it in one portion of the sample and not another would be high.

Site	FS col/100 mL	<i>C. perfringens</i> P/A	Salmonella (BSA) PEA	Salmonella (HEA) P/A
River water	63	yes	Absence	Presence
Creek Water	710000	yes	Presence	Absence
Rainwater	5	no	Absence	Absence

In summary, the best correlation of H₂S colour development time with other bacteria levels was for faecal coliforms, faecal streptococci and Clostridium perfringens. This is similar to other studies (see WHO 2002) and indicates the suitability of the H₂S test for testing drinking water for faecal pollution.

Field testing of the Hydrogen Sulphide (H₂S) test following Cyclone Ami

The difficulties in conducting water quality monitoring on remote islands are increased following natural disasters such as cyclones. An evaluation of the H₂S test against the conventional indicator organisms (total and faecal coliforms) was undertaken following the occurrence of Cyclone Ami in the Fiji Islands (SOPAC technical report #?). The H₂S test turned positive (black) for most samples that had faecal and total coliform levels above the respective WHO guidelines. Only about 11% and 8% of the samples that showed faecal and total coliform bacteria respectively did not test positive in the H₂S test, yielding 'false-negative' results. Similar disparities have been observed in other studies (see WHO 2002a for a summary) and are not unexpected as the coliform tests measure different bacterial groups than the H₂S test. Also the sample volume used in the H₂S test (10mL) is less than for the coliform-type indicators (100 mL), so the statistical probabilities of finding bacteria will differ. Lastly, there is an increased risk of introducing bacterial contamination when collecting and examining samples for faecal and total coliforms, due to the increased number of handling and filtration procedures, as compared to the H₂S test. This is particularly relevant when using the membrane filtration method in difficult non-laboratory conditions prevalent on outer islands. Upon closer examination of our results we found many of the 'false negative' samples had quite low levels of faecal and total coliform bacteria (e.g. see Labasa water depot sample results in Table 1 of SOPAC technical report #?). Therefore we believe that in many cases people drinking water that gave 'false negative' results in the H₂S tests would not necessarily be exposed to an increased risk to waterborne diseases.

About 2% and 6% percent of the samples that tested positive in the H₂S test did not have any total and faecal coliform bacteria respectively present. These are termed possible 'false-positives.' Similar findings have been documented in the literature (see WHO 2002 for a summary). This is likely due to the fact that some H₂S reducing bacteria (e.g. *Clostridium* sp.) persist in the environment longer than coliform bacteria. (WHO 2002). It could possibly be due to naturally occurring sulphide reducing bacteria being present, but the conditions needed for these bacteria to thrive are anaerobic waters with high organic matter and sulphate content. None of the waters we sampled fitted this description so we consider these results are unlikely to be false-positives in the sense of a natural H₂S producer being present. In any case, a false positive result indicates a problem, which when used in conjunction with a sanitary survey can provide information that would result in the suspect water either not being used, justify the system being cleaned and disinfected, the supply being disinfected or would suggest additional testing using conventional means. Nevertheless, following a disaster event, any positive test results should be regarded as unsafe for drinking purposes and disinfected.

We conducted a second visit to Vanua Levu, and several of the positive H₂S tests were re-tested for the presence of the spore-forming anaerobic bacteria, *Clostridium perfringens*, which is a strong H₂S producer and an indicator of faecal pollution. The results showed a relationship with the speed of H₂S development. Nearly 50% of the samples returned positive *C. perfringens* results, which indicated faecal contamination had entered into water supplies and/or the reticulation system. About 25% of samples that tested positive with the H₂S test had undetectable levels of *C. perfringens*. However, these were generally the samples where the H₂S test was slow to turn black, indicating that few H₂S producing micro-organisms were present when the sample was collected. Although it is suggested that further research on the quality of drinking waters in the Pacific should be performed to confirm the link between positive H₂S test results and the presence of faecal contamination and pathogens, the results thus far indicate that this test is valid.

Conclusion

The test is well suited for testing drinking water supplies for faecal contamination in the Pacific Islands, particularly in remote rural and outer island areas. The significant advantages of the H₂S test compared to other conventional microbiological (faecal and total coliforms) analyses is that it is very low in cost and does not require sophisticated equipment to manufacture or carry out the analyses. H₂S kits can easily be produced in Pacific Island countries where laboratories are often poorly equipped, or distributed by a regional organization, when needed. The results from H₂S tests are visual and therefore simple for people to understand, as a black colour change occurs when bacteria levels in drinking water are high. This enables communities and community health workers with minimum training to safely test their own water supplies. The time the H₂S test takes to turn black shows a correlation with faecal levels so an indication of the risk that pathogenic organisms are present can be obtained.

References

- APHA (1995). Standard Methods for the Examination of Water and Wastewater (1995). eds. Eaton, A., Clesceri, L., and Greenberg, A. 19th Edition. Washington, D.C. American Public Health Association, pp 9-19.
- Bartram, J. & Balance, R. eds. (1996). Water Quality Monitoring: A Practical Guide to the Design and Implementaiton of Freshwater Quality Studies and Monitoring Programmes. London: UNEP/WHO, Appendix 1.
- Castillo G, Duarte R, Ruiz Z, Marucic MT, Honorato B, Mercado R, et al. 1994. Evaluation of disinfected and treated drinking water supplies in Chile by the H₂S paper strip test. *Water Res* 28:1765-1770.
- Dillion, P. (1997). Groundwater pollution by sanitation on tropical islands. International Hydrological Programme, technical documents in hydrology no. 6. UNESCO, Paris.
- Dutka BJ. 1990. Review of the IDRC project to evaluate four simple, inexpensive, microbial water quality testing procedures. In: The use of simple, inexpensive microbial water quality tests: Results of a three-continent, eight-country research project (Dutka BJ, El-Shaarawi AH, eds). Rivers Research Branch, National Water Research Institute, Canada Centre for Inland Waters publication.
- Falkland T. 1999. Water resources issues of small island developing states. *Nat Resour Forum* 23:245-260.
- Manja KS, Maurya MS, Rao KM. 1982. A simple field test for the detection of faecal pollution in drinking water. *Bulletin World Health Organisation* 60:797-801.
- Martins M, Castillo G, Dutka BJ. 1997. Evaluation of drinking water treatment plant efficiency in microorganism removal by the coliphage, total coliform and H₂S paper strip tests. *Water Sci Technol* 35:403-407.
- Prüss, A., Kay, D., Fewtrell, L., & Bartram, J. (2002) Estimating the Burden of Disease from Water, Sanitation, and Hygiene at a Global Level. *Environmental Health Perspectives*. 110:537-542 (2002)
- WHO (1996). Guidelines for Drinking Water Quality, Second Edition, Volume 2. Geneva: World Health Organization, pp 82-88.
- WHO (2002). Evaluation of the H₂S for detection of faecal contamination in drinking water. WHO/SDE/WSH/02.08. World Health Organization, Geneva.

Appendix 1

Household Name:.....Phone:
 Street address/location:.....

Water Quality Testing Using the Hydrogen-Sulphide (paper strip) Test

Good water is essential for health. You can use this simple paper strip test to determine if your water is safe to drink. If the tube turns black within three days of collecting the sample, you must treat your water before drinking it, as harmful bacteria may be present.



When collecting a sample:

1. Allow the tap to run for about 30 seconds.
2. Uncap the lid of the sample tube and fill with water up to the mark.
3. Tightly cap the sample tube and put in a dark place to incubate for 3 days.
4. For comparison purposes, collect a sample of water that you know is safe, like boiled water or commercially bottled water.
5. Observe the samples every 12 hours (morning and evening) for 3 days and note below if water and /or the paper strip turns black. The water in the tube will have a slight brownish colour immediately after collecting the sample. This is due to the chemicals used and is considered normal.

DAY - 1		DAY - 2		DAY - 3	
TIME IN TUBE	12 hours	24 hours	36 hours	48 hours	60 hours
COLOUR CHANGE NOTED (Y/N)	<i>Yes/ No</i>	Yes/ No	Yes/ No	Yes/ No	Yes/ No
<i>High Risk</i>		<i>Moderate Risk</i>		<i>Low Risk</i>	

Note: If the sample tube turns black within 3 days of collection this indicated the presence of bacteria that could cause disease. In such cases the water should be treated before drinking. Treatment consists of boiling, filtering⁵, adding chlorine (household bleach), or exposing a 600ml clear plastic bottle of water to direct sunlight for a period of 4-6 hours.

Boiling means bringing the water to a rolling boil for a minimum period of 2 minutes.

Chlorine can be administered by adding 4-5 drops of household bleach to each litre of water treated or by adding chlorine tablets (meant for purifying drinking water) according to the instructions on the box, usually one tablet per litre.

⁵Filtering does not kill bacteria, it only reduces the number and lowers the risk to health.

For further information contact: SOPAC, Ph (679) 338 1377 or watersector@sopac.org

Appendix 2

WATER QUALITY TESTING - DATA SHEET

Address: _____

Location of source (describe): _____

SAMPLE NUMBER	TYPE WATER SOURCE: (Deep well - borehole; shallow well; surface water; spring; etc.)	DATE: of sample collection	TIME: of sample collection	LOCATION: (place where sample is collected)	RESULTS ² :							
					DAY 1		DAY 2		DAY 3			
					Date:	TIME	Date:	TIME	Date:	TIME		
				REMARKS ¹								
1												
2												
3												
4												
5												
6												
7												
8												
9												
10												
11												
12												

Notes: _____

1. Indicate under "remarks" if the water is visible turbid, colored, or contains settleable solids or material in suspension. Also, note any problem at the sampling site like a leaking tap, area unclean or littered, drainage problems, etc.
2. Results: a (-) indicates a negative; a (+), grey color, the reaction has started; (++) the reagent pad is now partially black; (+++) the reagent pad and the water is noticeably black.
3. Notes: Indicate the distance between the water source and any sources of pollution, like a compost pit, septic tank, leach field, etc.

Appendix 3: Rainwater Cistern Systems - Sanitary Survey

Country: _____ Date of visit: _____

Atoll: _____ Island: _____

Name of community/ village: _____ Est. pop. of village: _____ Est. # (households) in village: _____

Name of family (*it is suggested that the tank be numbered for easy identification*): _____

Number of people in family that drink water from this tank: _____ *[Note location of tank on map – use same number as sample]*

Sample No.: _____ *[Number that corresponds to number on H₂S sample tube]*

- 1. Condition of ROOF: good ⁽⁰⁾ _____; fair ⁽¹⁾ _____; poor ⁽²⁾ _____
- 2. Condition of guttering: good ⁽⁰⁾ _____; fair ⁽¹⁾ _____; poor ⁽²⁾ _____
- 3. Guttering sloped to drain: yes ⁽⁰⁾ _____; no ⁽²⁾ _____
- 4. Inlet screened or protected: yes ⁽⁰⁾ _____; no ⁽²⁾ _____
- 5. Interior tank clean: good ⁽⁰⁾ _____; fair ⁽¹⁾ _____; poor ⁽²⁾ _____
- 6. Condition of tank: good ⁽⁰⁾ _____; fair ⁽¹⁾ _____; poor ⁽²⁾ _____
- 7. Method of withdrawal by tap: yes ⁽⁰⁾ _____; no ⁽²⁾ _____
- 8. Tap and other plumbing in good repair: yes ⁽⁰⁾ _____; no ⁽²⁾ _____
- 9. Method for diverting first flush available: yes ⁽⁰⁾ _____; no ⁽²⁾ _____
- 10. No vegetation overhanging roof catchment area: yes ⁽⁰⁾ _____; no ⁽²⁾ _____

SCORE

TOTAL:

OTHER INFORMATION:

- 1. Type GUTTERING: vinyl _____; PVC _____; metal _____; other [describe]: _____

- 2. Type TANK: fibreglass _____; poured concrete _____; ferrocement _____; PVC _____; galvanized iron _____; other: [describe] _____

- 3. Location of tank: On raised platform _____; at ground level _____; partially below ground _____; majority of tank below ground _____
- 4. Estimated capacity of tank, in m³: _____ ($V=\pi r^2 h$)
- 5. Date constructed: _____ Date when last cleaned: _____
- 6. Other method of withdrawal [example, a bucket]: _____
- 7. Describe method for diverting first flush, if available (sketch on reverse): _____

RECOMMENDATIONS: _____

1. H₂S test (score): negative ⁽⁰⁾ _____; positive, day 1 ⁽¹¹⁾ _____; day 2 ⁽⁹⁾ _____; day 3 ⁽³⁾ _____

2. Sanitary Survey- Risk Score: _____

Relative risk = 1+2: LOW (<5) _____; MODERATE (5-10) _____; HIGH (>10) _____

Appendix 4: Cost of H₂S media

Recent prices in New Zealand dollars (quoted by Biolab Scientific) for the H₂S media

	Cost (NZD)	Cost per gram (NZD)
Peptone 500g	\$174.80	0.35
Di potassium hydrogen orthophosphate 500g	\$103.60	0.21
Ferric ammonium citrate 100g	\$30.20	0.30
Sodium Thiosulphate	\$56.99	

According to the formula;

40g peptone = \$14.00

3g Di potassium hydrogen orthophosphate = \$0.63

1.5g Ferric ammonium citrate = \$0.45

2g Sodium Thiosulphate = \$0.22

The total is: \$15.30 for enough chemicals to make 200 tests or roughly NZD \$0.08 per test.

This does not count the absorbent paper, sample container or lab time and equipment.

We can assume these could be free of charge. We obtain free pre-form bottles from Coca Cola company.

Compare this to the cost of an ampoule of MF-endo or other media, the H₂S test is much cheaper. Also, with the other tests you have real lab time and equipment use involved.

The H₂S can adequately indicate when a risk to human health occurs there is no excuse for not undertaking water testing because of a lack of financial or material resources.

As noted above: the test or chemicals may also be obtained on a 'cost of manufacture only' basis from the Institute of Applied Sciences, University of the South Pacific.



LIVE & LEARN ENVIRONMENTAL EDUCATION



World Health Organization



Australian Government
AusAID

